



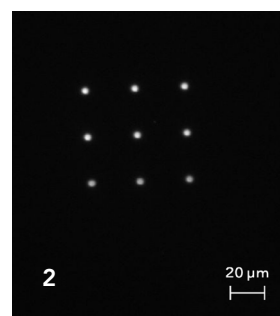

SPOTTING

FluidFM® spotting enables you to reproducibly create high density arrays with femtoliter spot sizes, both in gaseous and liquid environments. Print oligonucleotides, proteins, DNA, bacterial clones, oil and other materials with unprecedented precision and within their native environment. Applications of **FluidFM®** spotting are found in life sciences, sensor technology, minimal lubrication and beyond.

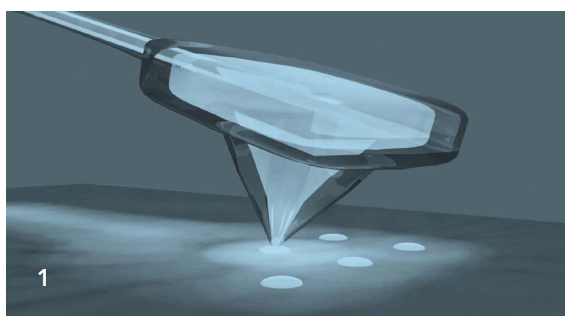

SPOTTING. 1 Illustration of **FluidFM®** nanopipette spotting. 2 Fluorescent image of femtoliter spots printed with **FluidFM®**. Courtesy of Behr P., ETH Zurich

PRECISE

fL SPOT DISPENSING

VERSATILE

LIQUID & GAS ENVIRONMENT

FAST

EXCHANGE OF INKS

LIMITLESS

CHOICE OF INKS

FluidFM® GIVES YOU THE EDGE.

FluidFM® based spotting redefines volumetric control of material spot printing. With **FluidFM®** spotting it is possible to reliably produce spot sizes with volumes as low as a few femtoliters on a large variety of target surfaces. The entire procedure can thereby be carried out in gaseous or liquid environment, with the option of additional humidity control. Precise adjustment of the relevant process parameters enables the creation of various spot sizes as required by your application.

Create microarrays of numerous biomolecules directly within their native environment to avoid denaturation of

your sensitive samples. Create functionalized surfaces with unparalleled spatial resolution for the creation of novel biosensing devices with smaller footprints and higher levels of performance. Provide minimally lubricated surface modifications for your most demanding micromechanical components. **FluidFM®** based spotting gives you the edge for your small scale material dispensing applications.

THE PROCEDURE IN BRIEF.

Spots are created while the **FluidFM®** nanopipette is in contact with the target surface via application of a short pressure pulse generated by the

FluidFM® microfluidics control system. Interaction forces with the surface are thereby monitored and adjusted in real-time during the entire procedure. Reproducible control of spot size and dispensing volumes can be comfortably achieved via modification of the contact time and applied pressure.

PUBLICATIONS

- 2009. A. Meister, J. Polesel-Maris, P. Niedermann, J. Przybylska, P. Studer, M. Gabi, P. Behr, T. Zambelli, M. Liley, J. Vörös & H. Heinzelmann. Nanoscale dispensing in liquid environment of streptavidin on a biotin-functionalized surface using hollow atomic force microscopy probes. *Microelectronic Engineering*, 86(4-6), 1481–1484. doi:10.1016/j.mee.2008.10.025
- 2014. H. Dermutz, R.R. Grüter, A.M. Truong, L. Demkó, J. Vörös & T. Zambelli. Local polymer replacement for neuron patterning and in situ neurite guidance. *Langmuir : the ACS journal of surfaces and colloids*, 30(23), 7037–46. doi:10.1021/la5012692

CONTACT US.

We offer complete support for our customers and distributors. Please visit the Cytosurge Help Center in order to access the **FluidFM®** user community. www.fluidfm.com

CYTOSURGE®

CYTOSURGE AG, SÄGEREISTRASSE 25, 8152 GLATTBRUGG, SWITZERLAND
PHONE +41 44 533 14 50, FAX +41 44 533 14 59, WWW.CYTOSURGE.COM

